AutoRep™ E Repeating Dispenser

Operating Manual

Please follow step by step!

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RAININ INSTRUMENT, LLC.

E-Mail: pipets@rainin.com

Internet: www.rainin.com

9978 24 · Printed in Germany · 13/1101

Mode d'emploi

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RAININ
®

Contents

	Page
Safety considerations	4
Purpose and limitations	7
Operating exclusions	8
Operating elements	9
Initial operation	10
Insert battery pack	10
Charge battery pack	11
Replace battery pack	14
Switching on/off	15
Inserting tips	16
Removing tips	18
Functions of the AutoRep E	19
Aspirating sample	20
DISP mode	22
AUTO DISP mode	24
PIP mode	26
Adjusting STEP volume	27
Adjusting aspiration and dispensing speed	28
Changing mode	30
Use of compatible tips from other manufacturers	31
Residual stroke	33
Non-contact dispensing	33
Maintenance and cleaning	33
Default settings	33
Troubleshooting	36
Technical data	42
Accuracy table	43
Ordering data	44
Return for repair	46

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Safety considerations

This instrument may sometimes be used with hazardous materials, operations, and equipment. To address all of the safety problems associated with its use in such applications is beyond the scope of this manual. It is the responsibility of the user of this instrument to consult and establish appropriate safety and health practices and determine the applicability of regulatory limitations prior to use.

1. Observe general instructions for hazard prevention and safety instructions; e.g. wear protective clothing, goggles and gloves. Observe standard laboratory procedures and precautions when pipetting infectious or hazardous samples.
2. Read and follow this Operating Manual.
3. Observe all safety precautions provided by reagent manufacturers. If in doubt, contact the manufacturer or supplier.
4. Use only with samples that do not attack polypropylene and polyethylene.

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Safety considerations

- continued -

5. Never use force on the instrument.
6. Operate only with tip attached (use Encode tips).
7. Avoid touching the tip orifice when working with hazardous samples.
8. Always work in a way which neither endangers the user nor any other person. Dispense samples into a suitable container, avoiding splashing.
9. Only press the STEP key after ensuring that dispensing the sample will not result in any danger. Dispense into a suitable vessel.
10. When a tip is inserted and locked the plunger can move inside the tip. If the tip was previously used, make sure that no sample was left inside, or that the tip is directed to a receiving vessel.
11. Do not use the instrument in potentially explosive environments or with explosive samples.
12. Only use the AC adapter indoors. Protect from humidity.

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Safety considerations

- continued -

13. To charge the NiMH batteries use only the original AC adapter in combination with the original charging dock.
14. Do not replace the original battery pack with other battery packs or cells. Use only original manufacturer's replacement battery packs.
15. Only dispose of completely discharged battery packs. Observe relevant disposal regulations for your community.
16. Use only original manufacturer's accessories and spare parts.
17. In case of malfunction, immediately stop dispensing. Consult the "Troubleshooting" section and call 800-543-4030 (in U.S.) if needed.

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Warning!

Improper use of the instrument or the batteries (short circuit, mechanical damage, overheating, incorrect AC adapter etc.) may, in extreme cases, lead to explosion of the batteries.

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Purpose and limitations

The AutoRep E is a microprocessor-controlled battery-operated repetitive pipette. The instrument automatically recognizes new Encode™ Syringe Tips, which are encoded for tip volume. This allows fastest possible volume selection. If you want to use compatible tips from other manufacturer, you must set the volume manually.

From now on, the term "tips" refers to Encode Syringe Tips and compatible tips.

The instrument is designed for repetitive dispensing of samples, observing the following limitations

- +15 °C to +40 °C (59 °F to 104 °F)
(instrument and reagents
- other temperatures on request)
- vapor pressure up to 500 mbar
- viscosity:
20 mPa s with 50 mL Encode tips
260 mPa s with 5 mL Encode tips
977 mPa s with 1.25 mL Encode tips

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Operating exclusions

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During proper operation, the dispensed sample only comes in contact with the tip and not with the AutoRep E. Never dispense samples which attack polypropylene or polyethylene.

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Do not use aggressive cleaning agents to clean the pipette.

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Note:

The instrument and the Encode tips are not autoclavable (121 °C / 250 °F). See sterile Encode tips on page 45.

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Initial operation

Insert battery pack

1. Open the battery compartment cover and insert battery pack, ensuring that the electrical contacts on the battery pack connect with those of the AutoRep E.
2. Replace the battery compartment cover.

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Warning!

The instrument must not be used or charged in potentially explosive environments.

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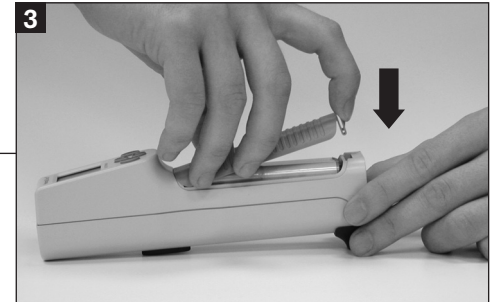
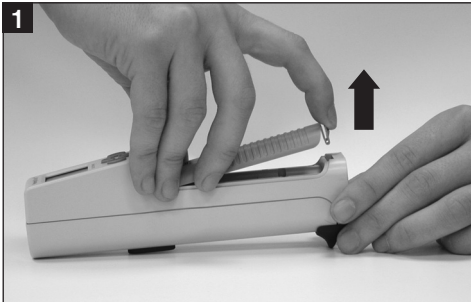
Charge battery pack

The battery pack can be charged either inside the AutoRep E or in the charging dock.

Before first use charge the battery pack inside the AutoRep E in the original charging dock for at least 3-½ hours. After 2-3 further charging cycles inside the instrument, the full capacity is reached.

Charge indicator is most accurate when battery is charged 3-½ hours in the instrument.

Removal of battery resets indicator. Upon reinsertion battery-charge frame flashes to indicate that charge status is approximate

Note:

Stored data and settings are retained when the battery pack is completely discharged or during replacement.

Remove the battery pack when not in regular or daily use.

- A tone indicates that the battery pack has to be charged or replaced. The battery symbol will flash on the display.

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Cargar dentro del AutoRep™ E

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Attention:

Before charging, make sure that the AC adapter is appropriate for your line voltage.

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1. Plug the AC adapter connector into the side of the charging dock.
2. Place the instrument with the battery pack into the charging dock. A tone indicates proper positioning of the instrument.
3. A green light on the charging dock indicates that the instrument is charging.

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The charging time for a nearly empty battery pack mounted in the instrument is approximately 3-½ hours.

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Attention:

Upon each battery insertion, charge-indicator frame flashes until battery is charged for at least 3-½ hours inside instrument.

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Charging the battery pack in the charging dock

The charging dock allows you to charge a separate battery pack while the AutoRep E is in use.

1. Plug the AC adapter connector into the side of the charging dock.
2. Insert the battery pack into the central opening of the charging dock, making sure the battery contacts connect with the charging dock contacts.
3. A green light on the charging dock indicates that the battery pack is charging.

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The charging time for a nearly empty battery pack in the charging dock is approximately 2-½ hours.

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Replace battery pack

A fully charged battery pack provides enough power for approximately 2-½ hours of continuous use with a 12.5 mL Encode Tip filled with a sample with similar viscosity and density to water.

- A tone indicates the need for recharging.
- The battery symbol will flash on the display.

When indicated, the battery should then be recharged or replaced. Before removing the battery, make sure the tip is completely empty.

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Switching on/off

- The AutoRep E can be switched on by inserting and locking a tip, or by pressing the ENTER key.
- When the locking lever is in the open position, the display shows "CLO" (Close).
- Display: **rEF** (reference run): Press the STEP-key to start reference run.
rEF is displayed after every replacement of the battery pack.

- Auto Power Off

- The AutoRep E is factory-set to switch off after 3 minutes of inactivity.
- To switch off manually, press the ENTER key briefly

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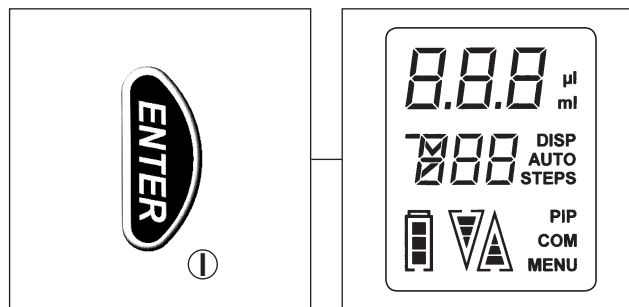
Inserting tips

Attention:

Risk of injury! Do not touch the plunger seat inside the instrument!

1. Push the locking lever to its upper stop.
2. Slide a tip laterally into the AutoRep E.
3. Push the locking lever down until you hear a distinct clicking (2 steps).
4. If using Encode tips, the tip capacity is recognized by the AutoRep E and its capacity is shown briefly on the display.
5. Then the partial volume to be dispensed in each step is displayed. The default setting is 10 % of the maximum tip capacity.

If using compatible tips (not Encode tips) you must enter the capacity of the tip yourself.

Note:

When using 25 mL and 50 mL Encode Tip, the supplied adapter must be used. It can be connected to the Encode tips using a twist lock and can be removed after use.

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

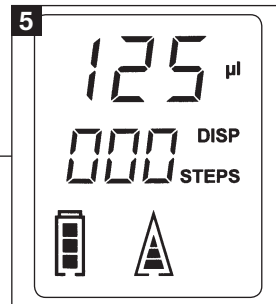
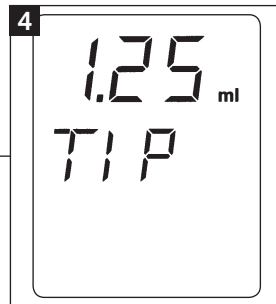
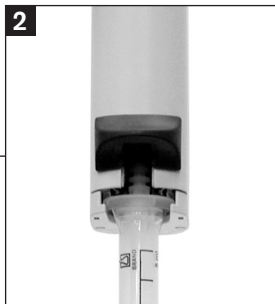
- If the locking lever is not fully engaged, horizontal bars are shown on the display.
- If the locking lever was fully closed but no tip was inserted, **"no TIP"** appears on the display.

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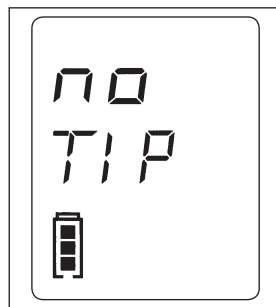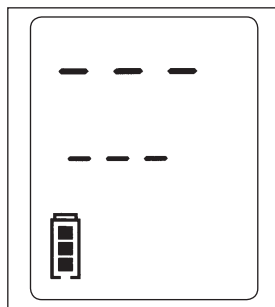Removing tips

1. Hold the AutoRep E over a suitable container.
2. Completely discharge the tip by continuously pressing the STEP key (releasing the residual stroke lock), or by using the "E" mode (see page 23).
3. Push the locking lever to its upper stop.
4. Remove the tip.

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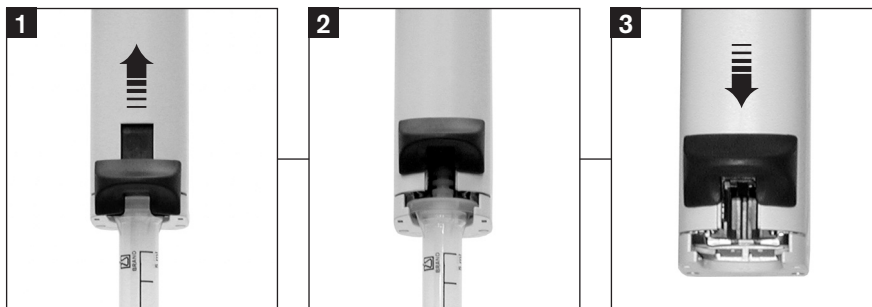
Note:

After dispensing highly viscous or crystallizing sample, the Encode tip may wear and require replacement

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Functions of the AutoRep™ E

The instrument performs the following functions:

1. Dispensing **DISP** (see page 22): Default mode after switching on. A user-defined volume can be dispensed repeatedly, depending on the capacity of the tip and the filling volume.
2. Automatic Dispensing **AUTO DISP** (see page 24): Identical to DISP mode, but with automatic dispensing by continuously pressing the STEP key. Releasing the STEP key interrupts the process.
In the AUTO DISP mode the instrument has a "**Learn**" function: After pipetting at least three steps, the AutoRep E calculates the average dispensing interval, and will repeatedly dispense at that interval as long as the STEP key is pressed.
3. Pipetting **PIP** (see page 26): In "PIP" mode, the AutoRep E functions as a positive displacement non-repeating pipette, particularly useful for dense, volatile or viscous materials. A preselected volume is aspirated and dispensed.

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1. Push the STEP key once. The tip will be filled completely.
2. The arrow on the display points upward (Δ "Aspirate").

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Note:

To interrupt the filling process, press the STEP key. To continue filling, press the Δ key first and hold down the STEP key in succession. The step count on the display will increase.

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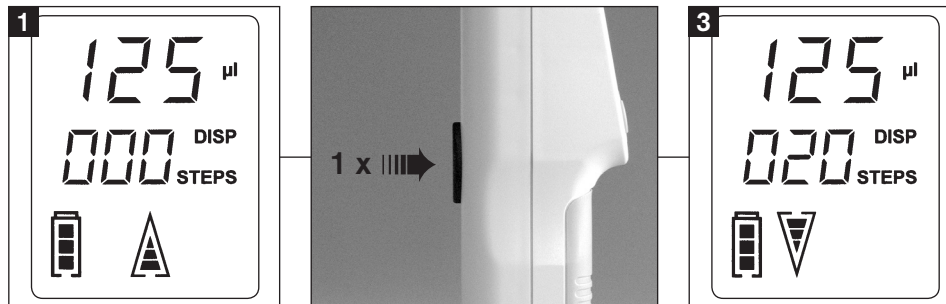
3. During filling, the number of remaining steps is continuously displayed.

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Attention:

After aspirating liquid, the Auto-Rep E will automatically dispense a small amount of liquid in order to compensate for gear backlash.

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DISP mode

1. The arrow on the display points downward (▽ "Dispense").
2. Press the STEP key to dispense the selected step volume (see "Adjusting STEP volume" page 27).
3. The display shows the remaining number of steps.
4. After dispensing the last full step, a small amount of sample will be left in the tip. To discharge this volume, hold down the STEP key.
5. Remove tip (see page 18).

Note:

The tip must be completely empty before you try to remove it.

Attention:

Dispensing can be interrupted at any time by pressing the Δ key.

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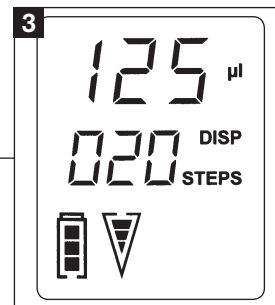
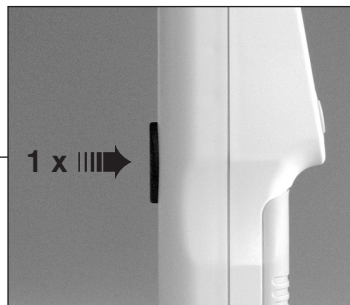
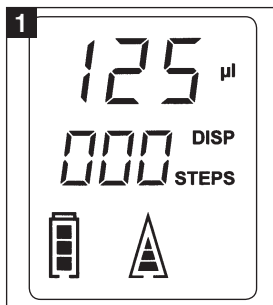
Note:

To empty the tip at any time, press the ▽ key and then hold down the STEP key. The letter E appears on the display. This process can be interrupted at any time by pressing the Δ key.

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A partially filled tip can be completely filled by pressing and holding, in sequence, the Δ key and the STEP key.

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⇒ siehe Seite 26.

AUTO DISP mode

1. To activate the AUTO DISP mode, briefly press the MENU key twice.

Note:

The AUTO DISP mode must be actively selected.

2. Select the AUTO DISP mode with the $\nabla\Delta$ keys (ON/OFF).
3. Confirm by pressing ENTER.
4. **"Learn"** feature: After pipetting at least three steps, the instrument automatically determines the average time interval between steps (see page 19).
5. Hold down the STEP key to dispense the selected volume (see page 27) at the calculated intervals.

Note:

Intervals can range from 0.1 to 5 seconds.

6. To interrupt dispensing, release the STEP key. The calculated step interval remains in memory.
7. In AUTO DISP mode, the "Learn" process can be restarted at any time.

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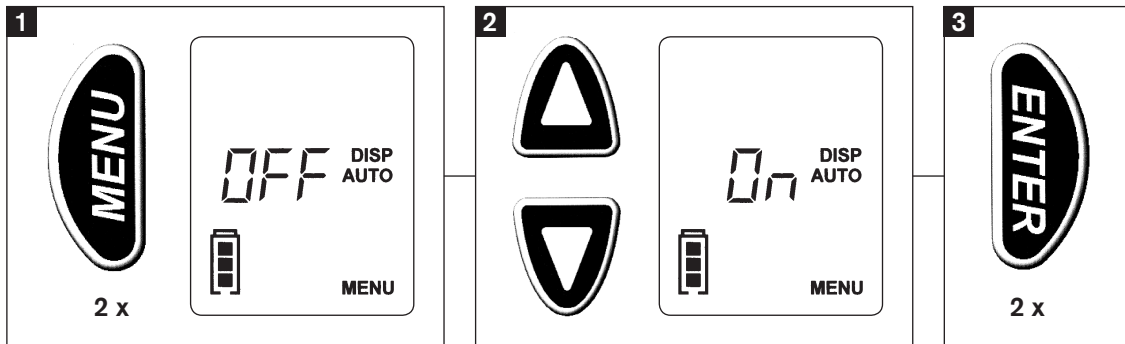
PIP mode

(Reverse pipetting, positive displacement principle)

1. Select the PIP mode by holding down the MENU key (> 3s), and press it briefly (see page 30).
Make sure that plunger is at lowest position!
2. Select a volume (page 27).
3. To aspirate the selected volume, press the STEP key. A volume slightly larger than the selected volume will be drawn into the tip.
4. To dispense the volume, press the STEP key again. Some residual volume will remain in the tip.
5. To aspirate a new sample, press the STEP key again.
6. When you have completed pipetting, discharge the residual stroke by holding down the STEP key.

Attention!

To interrupt the dispensing process at any time, press the Δ key.

Note:

To change from PIP mode to AUTO DISP mode, you have to change to DISP mode first.

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Adjusting STEP volume

Note:

After the Encode tip size has been recognized, the dispensing volume is shown on the display. The default step volume is 10 % of the nominal capacity.

1. Press the MENU key once. The volume (VOL) is shown on the display.
2. Use the $\nabla\Delta$ keys to select the appropriate volume. Holding down the $\nabla\Delta$ keys will accelerate the rate of change.
3. Confirm the volume selection by pressing ENTER.

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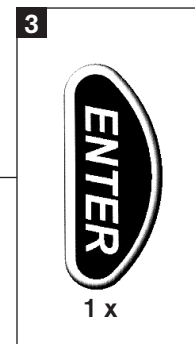
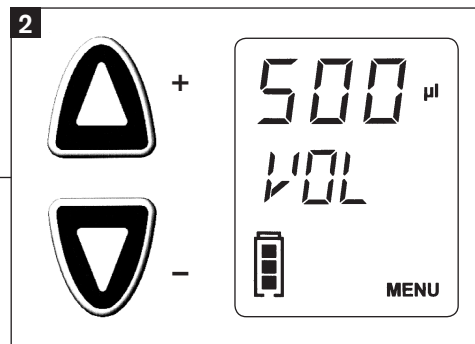
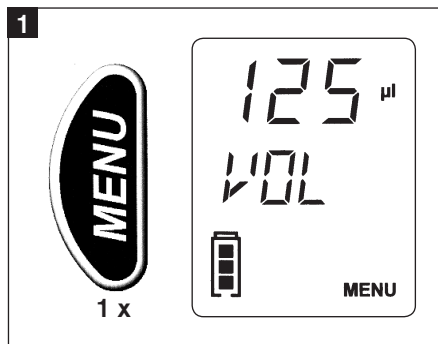
Adjusting aspiration and dispensing speed

1. Hold down the MENU key (> 3s) to adjust the aspiration speed. Choose between three different aspiration speeds.

Note:

Lowest speed:

Display shows outline of arrow with one bar.

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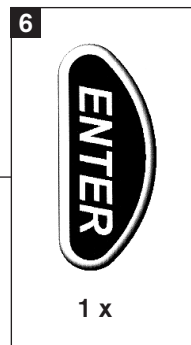
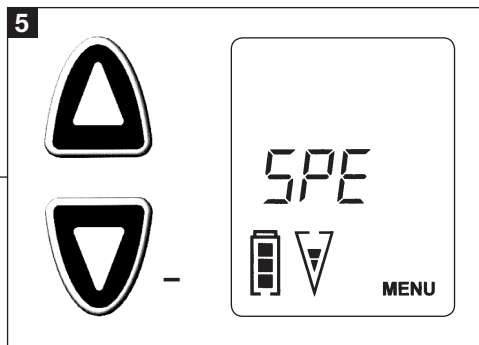
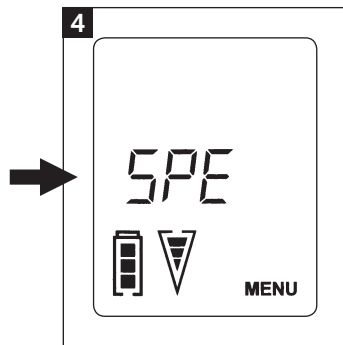
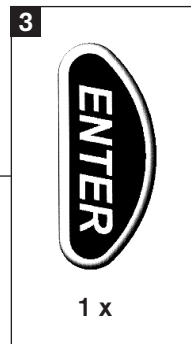
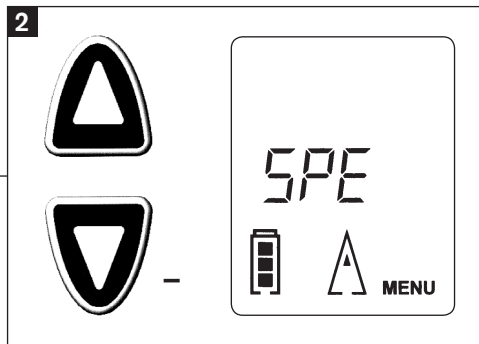
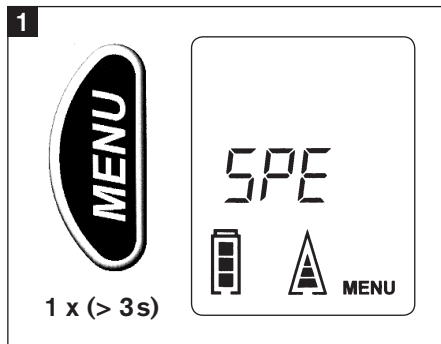
2. Use the $\nabla\Delta$ keys to select aspiration speed.
3. Confirm aspiration speed setting by pressing ENTER.
4. The down arrow will now flash.
5. Use the $\nabla\Delta$ keys to select dispensing speed.
6. Confirm settings by pressing ENTER. You are returned to the main menu.

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Changement de mode

Cambio de modo

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1. Hold down the MENU key (> 3s) and then press it again briefly. The mode can then be selected.

Make sure that plunger is at lowest position!

- Modes available are PIP and DISP.
- DISP is the default mode.

2. Use the $\nabla\Delta$ keys to select desired mode.
3. Confirm setting by pressing ENTER.

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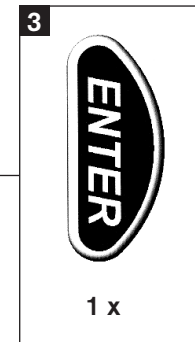
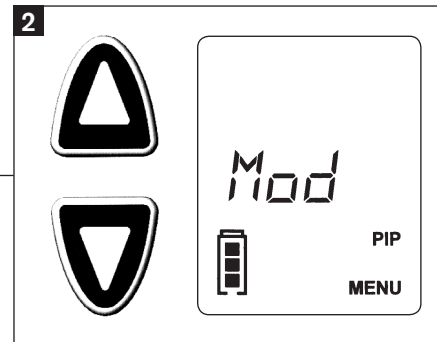
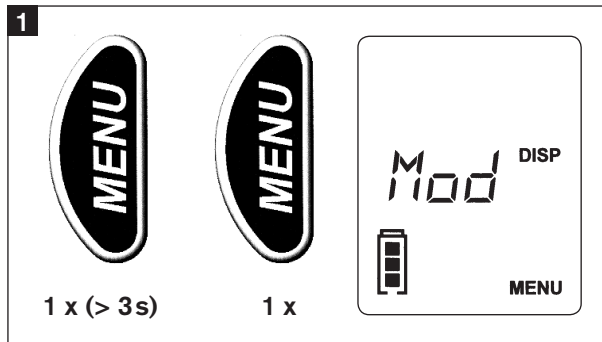
Use of compatible tips from other manufacturers

Compatible dispenser tips that are not encoded for size can be used with the AutoRep E (see "Insert tips" page 16). Bars are shown on the display to indicate an un-encoded tip. After inserting one of these tips, the display will automatically go to tip selection mode.

By default, the 5 mL tip is displayed flashing. 12 possible tip sizes are available.

1. Use the $\nabla\Delta$ keys to select appropriate tip size.
2. Confirm setting by pressing ENTER.

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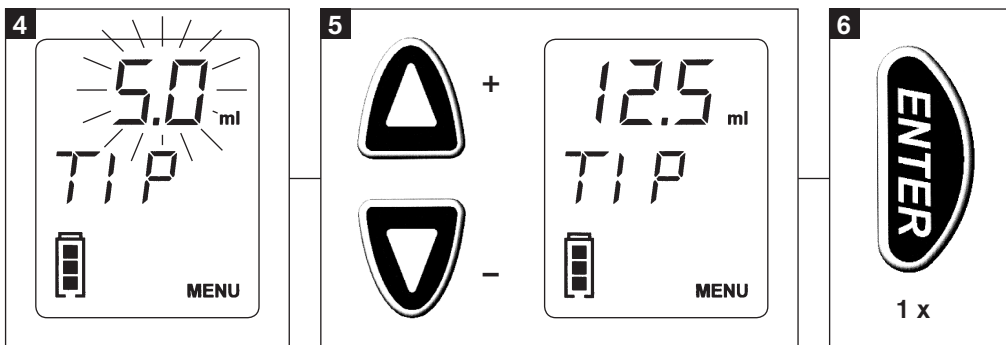
Compatible tips from other manufacturers

– continued –

Attention!

For convenience, when inserting the next unencoded tip, the previously selected tip size will come up first on the menu. Confirm tip size by pressing ENTER.

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At the time of printing, the following dispenser tips were compatible with the AutoRep E:

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Residual stroke

A lock prevents the residual stroke from being dispensed as a regular dispensed or pipetted volume. The symbol Δ indicates this remaining volume. It can be discharged by holding the STEP key.

Non-contact dispensing

When tips with volumes of 5 mL or larger are selected, the dispensing speed adjusts to allow accurate dispensing without contacting the walls of vessel (free-jet dispensing). This can be overridden by selecting a different speed (see p. 28).

Maintenance and cleaning

The AutoRep E is calibrated at the factory and is maintenance-free. If required, clean the instrument with a cloth dampened with water or a mild soap solution. The AutoRep E must not be dismantled.

Default settings

To revert the AutoRep E to its factory default settings, simultaneously press the MENU and ENTER keys for > 3 seconds, with the plunger in the lowest position.

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Troubleshooting

Problem	Error message displayed	Possible cause	Corrective action
Tip dripping	No message	Leaking tip	Replace tip
Faint display; reduced speed	Battery symbol flashing	Battery low	Recharge battery or replace with charged battery
Instrument not responding	ERR 001	<ul style="list-style-type: none"> - Locking lever not engaged during intake or dispensing - Locking lever open at low third of the tip 	<p>1. Press ENTER to confirm error.</p> <p>Follow instructions to avoid instrument damage!</p> <p>Display: CLO (Close)</p> <p>2. Engage the locking lever</p> <p>Display: rEF (Reference run)</p> <p>3. Press the STEP key</p> <p>Attention: Liquid remaining in tip will be discharged during the reference run (rEF). Discharge sample into a suitable vessel (see pages 4-6).</p>
Instrument not responding	ERR 002	<p>Motor blocked</p> <ul style="list-style-type: none"> - Battery low - Viscosity is too high (see page 7) - Tip orifice obstructed 	<p>1. Press ENTER to confirm error.</p> <p>Follow instructions to avoid instrument damage!</p> <ul style="list-style-type: none"> - Recharge or replace battery - Reduce intake/dispensing speed <p>Display: rEF (Reference run)</p> <p>2. Press the STEP key</p> <p>Attention: Liquid remaining in tip will be discharged during the reference run (rEF). Discharge sample into a suitable vessel (see pages 4-6).</p>
Instrument not responding	ERR TIP	Locking lever open at lower part of the tip	<p>1. Press ENTER to confirm error.</p> <p>Follow instructions to avoid instrument damage!</p> <p>2. Display: OPE (Open)</p> <p>Push the locking lever to its upper stop.</p>

Troubleshooting - continued -

Problem	Error message displayed	Possible cause	Corrective action
			3. Display alternating between out TIP and CLO Remove tip and close locking lever (without tip). 4. AutoRep E automatically carries out reference run. 5. Display: no TIP Insert new tip (see page 16)
Instrument not responding	ERR 003	Electronic fault	Press ENTER to confirm error. Send instrument in for repair.
No tone, no light when placing instrument into the charging dock	ACC	- Instrument contains no battery - Battery inserted improperly - Battery faulty	- Insert battery pack. - Make sure battery contacts are aligned with contacts in the instrument - Replace battery
Charge-indicator is flashing	No message	Removal or insertion of a new battery	- Charge battery for at least 3-½ hours inside the AutoRep E

Note:

If other error messages are displayed, remove battery pack and insert it again, or send instrument in for repair.

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Technical data

Battery pack:

Rechargeable Nickel Metal Hydride battery pack:
4 NiMH, AAA cells,
4.8 V/700 mAh

Charging times for battery:

in battery compartment of charging dock: approx. 2-½ h
inside the AutoRep E: approx. 3-½ h
When battery capacity starts to diminish, recharge battery in the battery compartment of the charging dock.

Battery capacity:

A full charge gives approx. 2-½ hours of continuous work with a 12.5 mL Encode Tip.

This sign certifies that the product meets the requirements of the EEC directive and has been tested according to the specified test methods.

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Ref.: ENC-SET

Cat. No.: ENC-SET

Adapter for 25 and 50 mL Encode tips, polypropylene, autoclavable

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Zur Reparatur einsenden

Important:

For safety reasons, only clean and decontaminated instruments will be accepted for inspection or repair.

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- Thoroughly clean and decontaminate the instrument.
- Call 800-543-4030 for a Return Authorization Number
- Send the AutoRep E in a suitable packing container to Rainin Service, Rainin Road, Woburn, MA 01801, describing:
- The nature of the problem
- What samples have been used

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Note:

No warranty claims can be considered if non-authorized repairs (repairs other than by trained Rainin technicians) have been made or attempted.

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