repetman®

Gebrauchsanleitung
User's Guide
Mode d'emploi
Instrucciones de manejo

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Please read carefully before use, and follow all operating and safety instructions!

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TABLE OF CONTENTS

SAFETY INSTRUCTIONS 4
PURPOSE AND LIMITATIONS OF USE 7
OPERATING EXCLUSIONS 8
OPERATING ELEMENTS 9
INITIAL OPERATION 10
Insert the battery pack 10
Charge the battery pack 11
Charge REPETMAN 12
Replace the battery pack 14
SWITCHING ON/OFF 15
INSERTING TIPS 16
REMOVING TIPS 18
FUNCTIONS OF REPETMAN 19
ASPIRATING LIQUID 20
DISP MODE 22
AUTO-DISP MODE 24
PIP MODE 26
ADJUSTING STEP VOLUME 27
ADJUSTING ASPIRATION AND DISPENSING SPEED 28
CHANGING MODE 30
USE OF COMPATIBLE TIPS FROM OTHER MANUFACTURERS 31
RESIDUAL STROKE 33
FREE –JET DISPENSING 33
MAINTENANCE AND CLEANING 33
DEFAULT SETTINGS 33
TROUBLESHOOTING 36
TECHNICAL DATA 42
ACCURACY TABLE 43
ORDERING DATA 44
RETURN FOR REPAIR 46
DISPOSAL 47
WARRANTY 48

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SAFETY INSTRUCTIONS

This instrument may sometimes be used with hazardous materials, operations, and equipment. To address all of the safety problems associated with its use in such applications is beyond the scope of this manual. It is the responsibility of the user of this instrument to consult and establish appropriate safety and health practices and determine the applicability of regulatory limitations prior to use.

1. Every user must read and understand this operating manual prior to using the instrument and observe these instructions during use.
2. Observe general instructions for hazard prevention and safety instructions; e.g. wear protective clothing, goggles and gloves. When working with infectious or other hazardous samples, all appropriate regulations and precautions must be followed.
3. Observe all safety precautions provided by reagent manufacturers.
4. Never use the instrument in an atmosphere with danger of explosion. Highly flammable liquids must not be pipetted.

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SAFETY INSTRUCTIONS - CONTINUED -

- Use the instrument only for dispensing liquids, with strict regard to the defined Operating Exclusions and Limitations. Observe operating exclusions (refer to page 7 and 8). If in doubt, contact the manufacturer or supplier.
- Always work in a way which neither endangers the user nor any other person. Avoid splashes. Use only suitable containers.
- When a tip is inserted and locked the plunger can move inside the tip. If the tip was previously used, make sure that no sample was left inside, or that the tip is directed to a receiving vessel.
- Only press the **STEP** key after ensuring that dispensing the sample will not result in any danger.
- Avoid touching the tip orifice when working with hazardous samples.
- Never use force on the instrument.
- Use only original manufacturer's accessories and spare parts. Do not attempt to make any technical alterations. Do not dismantle the instrument.
- Always check the instrument for visible damage before use. If there is a sign of a potential malfunction, immediately stop dispensing. Consult the **TROUBLESHOOTING** section of this manual on page 36 and 37, and contact the manufacturer if needed.

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SAFETY INSTRUCTIONS

- CONTINUED -

- The original battery must not be replaced with non rechargeable batteries or rechargeable batteries of other manufacturers.
- To charge the NiMH battery pack use only the original AC adapter in combination with the original charging dock.
- The AC adapter has to be protected against moisture and must be used only for this instrument.
- Dispose of battery pack only when completely discharged, and according to applicable regulations.

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Note:
Attention! Risk of injury.
Do not touch the plunger seat inside
the instrument!

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Warning!
Improper use of the instrument
or the batteries (short circuit,
mechanical damage, overheating,
incorrect AC adapter etc.) may, in
extreme cases, lead to explosion of
the batteries.

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PURPOSE

REPETMAN® is a microprocessor controlled battery—operated repetitive pipette. The instrument automatically recognizes new Gilson REPET-TIPS, which are encoded for tip volume. This allows fastest possible volume selection. If you want to use compatible tips from other manufacturers, you must set the volume manually.

From now on, the term “tips” refers to REPET-TIPS with size encoding and compatible tips.

LIMITATIONS OF USE

The instrument is designed for quick and simple repetitive dispensing of samples, observing the following limitations:

- +15°C to +40°C (59°F to 104°F) (instrument and reagents, other temperatures on request).
- Vapor pressure up to 500 mbar.
- Viscosity:

20 mPa.s with 50 mL REPET-TIPS
260 mPa.s with 5 mL REPET-TIPS
977 mPa.s with 1.25 mL REPET-TIPS

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OPERATING EXCLUSIONS

During proper operation, the dispensed sample only comes in contact with the tip and not with REPETMAN.

The user has to ensure the compatibility of the instrument with the intended application.

Never dispense samples which attack polypropylene, polyethylene or polycarbonate (housing). Avoid aggressive fumes (danger of corrosion)!

Note:

The instrument and the tips are not autoclavable (121°C / 250°F). See Positive Displacement Tips, on page 45. Pipette tips are disposable. For disinfection you may use commercially available disinfecting solutions as long as they do not attack PC/PBT or damage the display.

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INITIAL OPERATION

Insert the Battery Pack

1. Open the battery compartment cover and insert battery pack, ensuring that the electrical contacts on the battery pack connect with those of the REPETMAN.
2. Replace the battery compartment cover.

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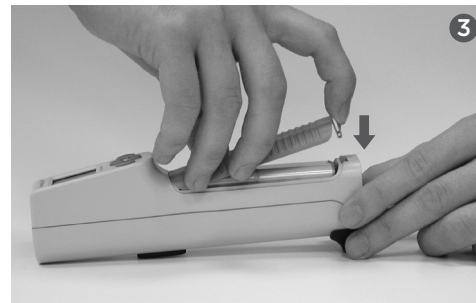
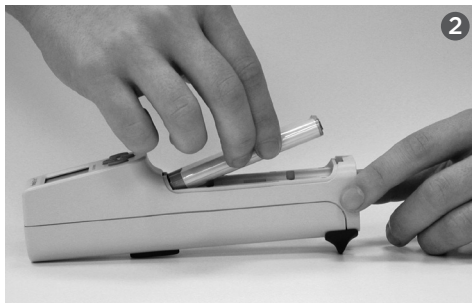
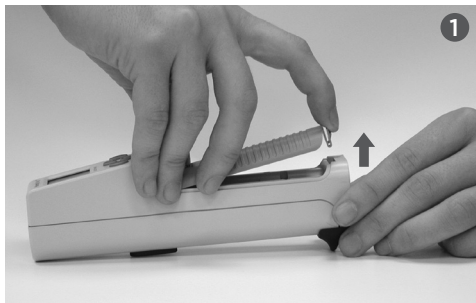
Warning!
The instrument must not be used or charged in potentially explosive environments.

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Charge the Battery Pack

The battery pack can be charged either inside the REPETMAN or in the charging dock.

Before first use charge the battery pack inside the REPETMAN in the original charging dock for at least 3.5 hours. After 2-3 further charging cycles inside the instrument, the full capacity is reached.

Charge indicator is most accurate when battery pack is charged 3.5 hours in the instrument.

Removal of battery pack resets indicator. Upon reinsertion battery charge frame flashes to indicate that charge status is approximate.

Note:

Stored data and settings are retained when the battery pack is completely discharged or during replacement. Remove the battery pack when not in regular or daily use.

- A single tone indicates that the battery pack has to be charged or replaced. The battery symbol will flash on the display.

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Charge REPETMAN

Caution!

Before charging, make sure that the AC adapter is appropriate for your line voltage.

1. Plug the AC adapter connector into the side of the charging dock.
2. Place the instrument with the battery pack into the charging dock. A tone indicates proper positioning of the instrument.
3. A green light on the charging dock indicates that the instrument is charging.

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The charging time for a nearly empty battery pack mounted in the instrument is approximately 3.5 hours.

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Caution!

Upon each battery pack insertion, charge—indicator frame flashes until battery pack is charged for at least 3.5 hours inside instrument.

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CHARGING THE BATTERY PACK IN THE CHARGING DOCK

The charging dock allows you to charge a separate battery pack while REPETMAN is in use.

1. Plug the AC adapter connector into the side of the charging dock.
2. Insert the battery pack into the central opening of the charging dock, making sure the battery pack contacts connect with the charging dock contacts.
3. A green light on the charging dock indicates that the battery pack is charging.

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The charging time for a nearly empty battery pack in the charging dock is approximately 2.5 hours.

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Replace the Battery Pack

A fully charged battery pack provides enough power for approximately 2.5 hours of continuous use with a 12.5 mL Repet-Tip filled with a sample with similar viscosity and density to water.

- A tone indicates the need for recharging.
 - The battery symbol will flash on the display.
- This indicates the need to recharge REPETMAN or to insert a spare battery pack.

Before removing the battery, make sure the tip is completely empty.

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SWITCHING ON/OFF

- REPETMAN can be switched on by inserting and locking a tip, or by pressing the **ENTER** key.
- When the locking lever is in the open position, the display shows “CLO” (Close).
- Display: **rEF** (reference run): Press the **STEP** key to start reference run. **rEF** is displayed after every replacement of the battery pack.

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Auto Power Off

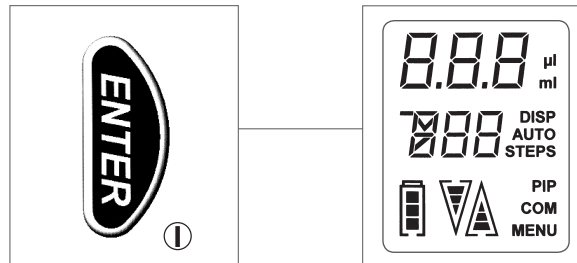
- REPETMAN is factory-set to switch off after 3 minutes of inactivity.
- To switch off manually, press the **ENTER** key briefly

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INSERTING THE TIPS

Caution!

Risk of injury! Do not touch the plunger seat inside the instrument!

1. Push the locking lever to its upper stop.
2. Slide a tip laterally into REPETMAN .
3. Push the locking lever down until you hear a distinct clicking (2 steps).
4. If using encoded REPET-TIPS from Gilson, the REPET-TIP capacity and the word **TIP** appear on the display for 3 sec.
5. Then the partial volume to be dispensed in each step is displayed. The default setting is 10% of the maximum tip capacity.

If using compatible tips (not REPET-TIPS) you must enter the capacity of the tip yourself (see page 31).

Note:

When using 25 mL and 50 mL tips, the supplied adapter must be used. It can be connected to the tips using a twist lock and can be removed after use.

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

6. If the locking lever is not fully engaged, horizontal bars are shown on the display.
7. If the locking lever was fully closed but no tip was inserted, **no tip** appears on the display.

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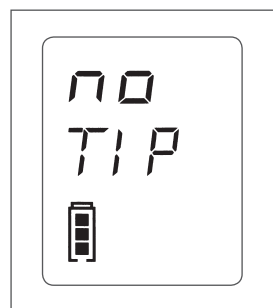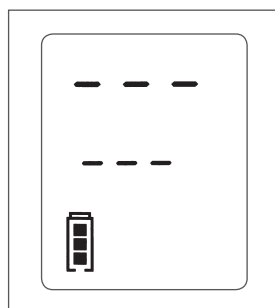REMOVING TIPS

1. Hold REPETMAN over a suitable container.
2. Completely discharge the tip by continuously pressing the **STEP** key (releasing the residual stroke lock), or by using the "E" mode (see page 23).
3. Push the locking lever to its upper stop.
4. Remove the tip.

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Caution!

Don't remove REPET-TIP before piston of the tip is in the lowest position (tip completely discharged!).

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Note:

After dispensing highly viscous or crystallizing sample, the tip may wear and require replacement.

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FUNCTIONS OF REPETMAN

The instrument performs the following functions:

1. Dispensing **DISP** (see page 22): Default mode after switching on. A user-defined volume can be dispensed repeatedly, depending on the capacity of the tip and the filling volume.
2. Automatic Dispensing **AUTO-DISP** (see page 24): Identical to **DISP** mode, but with automatic dispensing by continuously pressing the **STEP** key. Releasing the **STEP** key interrupts the process.
3. In the **AUTO-DISP** mode the instrument has a "Learn" function: After pipetting at least three steps, REPETMAN calculates the average dispensing interval, and will repeatedly dispense at that interval as long as the **STEP** key is pressed.
4. Pipetting **PIP** (see page 26): In **PIP** mode, REPETMAN functions as a positive displacement non-repeating pipette, particularly useful for volatile, foamy or high viscous materials. A preselected volume is aspirated and dispensed (Positive displacement principle).

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ASPIRATE THE LIQUID

Note:

Default setting for the REPETMAN is DISP mode (see page 22).

We recommend that a new REPET-TIP be bled before use. To this end, the tip should at a minimum be filled with the liquid and then emptied again (see page 21). Small air bubbles in the vicinity of the piston after bleeding do not affect the results.

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1. The arrow on the display points upward Δ “Aspirate”.
2. Push the **STEP** key once. The tip will be filled completely. During filling, the number of remaining steps is continuously displayed.
3. After the filling procedure has been completed, the arrow “Dispense” appears on the display.

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Note:

To interrupt the filling procedure and to bleed the **REPET-TIP**:

To interrupt the filling process, press the **STEP** key. To continue filling, press the Δ key first and hold down the **STEP** key in succession. The step count on the display will increase.

Activate the ∇ button and press and hold down the **STEP** button to completely discharge the liquid again.

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Caution!

After aspirating liquid, REPETMAN will automatically dispense a small amount of liquid in order to compensate for gear backlash.

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DISP MODE

1. The arrow on the display points downward ▽ “Dispense”.
2. Press the **STEP** key to dispense the selected step volume (see “Adjusting **STEP** volume” page 27).
3. The display shows the remaining number of steps.
4. After dispensing the last full step, a small amount of sample will be left in the tip. To discharge this volume, hold down the **STEP** key.
5. Remove tip (see page 18).

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Note:

The tip must be completely empty before you remove it.

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Caution!

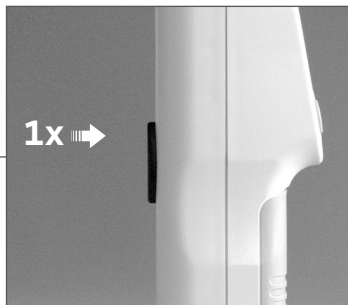
Dispensing can be interrupted at any time by pressing the Δ key.

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Note:

To empty the tip at any time, press the ▽ key and then hold down the STEP key. The letter E appears on the display and the arrow pointing downwards flashes. This process can be interrupted at any time by pressing the Δ key.

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A partially filled tip can be completely filled by pressing and holding, in sequence, the Δ key and the STEP key.

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AUTO-DISP mode:
• see page 24,
PIP mode:
• see page 26.

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AUTO-DISP-MODE

1. To activate the **AUTO-DISP** mode, briefly press the **MENU** key twice.

Note:

The **AUTO-DISP** mode must be manually selected.

2. Select the **AUTO-DISP** mode with the Δ keys (**ON/OFF**).
3. Confirm by pressing **ENTER**.
4. "Learn" feature: after pipetting at least three steps, REPETMAN automatically determines the average time interval between steps (see page 19).
5. Hold down the **STEP** key to dispense the selected volume (see page 27) at the calculated intervals.

Note:

Intervals can range from 0.1s to 5s.

6. To interrupt dispensing, release the **STEP** key. The calculated step interval remains in memory.
7. In **AUTO-DISP** mode, the "Learn" process can be restarted at any time.

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PIP MODE

(Reverse pipetting, positive displacement principle)

1. Select the **PIP** mode by holding down the **MENU** key (>3s), and press it briefly (page 30).
2. Make sure that plunger is at lowest position!
3. Select a volume (page 27).
4. To aspirate the selected volume, press the **STEP** key. A volume slightly larger than the selected volume will be drawn into the tip.
5. To dispense the volume, press the **STEP** key again. Some residual volume will remain in the tip.
6. To aspirate a new sample, press the **STEP** key again.
7. When you have completed pipetting, discharge the residual stroke by holding down the **STEP** key.

Caution!

To interrupt the pipetting process at any time, press the **Δ** key.

Note:

To change from **PIP** mode to **AUTO-DISP** mode, you have to switch to **DISP** mode first.

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ADJUSTING STEP VOLUME

Note:

After the tip size has been recognized, the dispensing volume is shown on the display. The default step volume is 10% of the nominal capacity.

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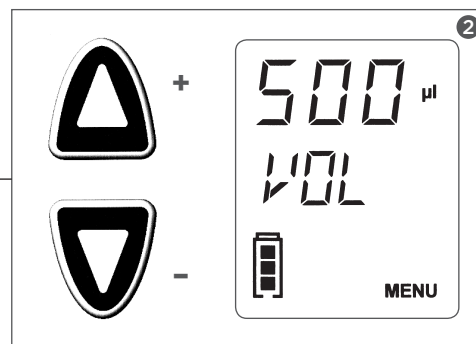
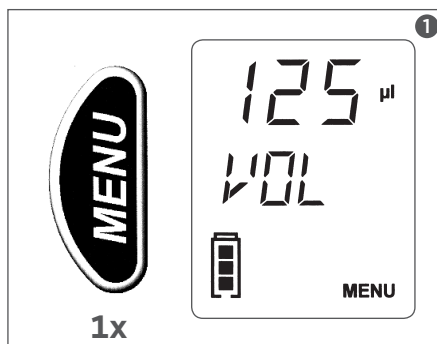
1. Press the **MENU** key once. The volume (**VOL**) is shown on the display.
2. Use the ∇ keys to select the appropriate volume. Holding down the ∇ keys will accelerate the rate of change.
3. Confirm the volume selection by pressing **ENTER**.

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ADJUSTING ASPIRATION AND DISPENSING SPEED

1. Hold down the **MENU** key (>3s) to adjust the aspiration speed. The arrow for selection of aspiration speed lights up on the display. Choose between three different aspiration speeds.

Note:

Lowest speed:
Display shows outline of arrow with one bar.

Highest speed:
Display shows outline of arrow with three bars.

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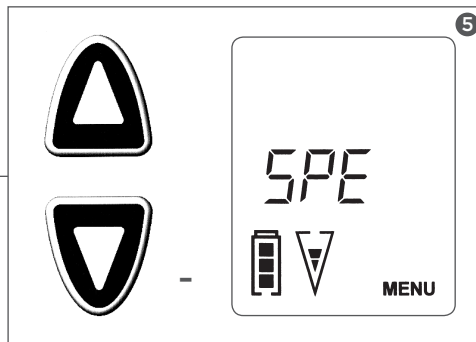
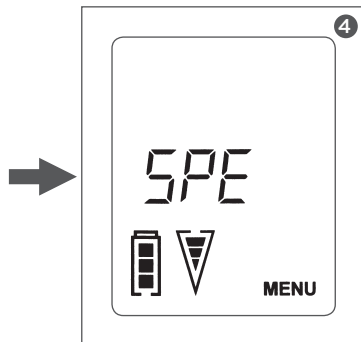
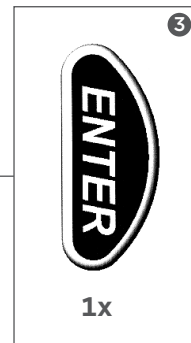
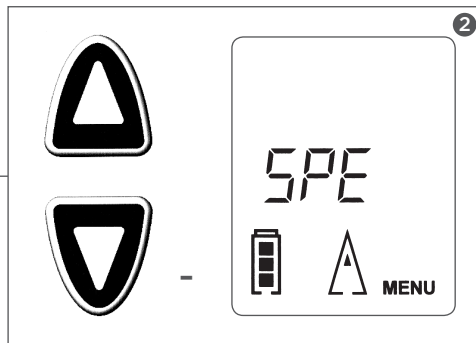
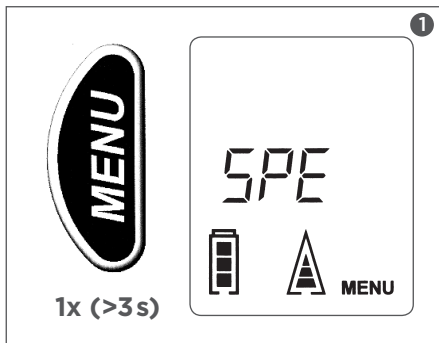
2. Use the ∇ keys to select aspiration speed.
3. Confirm aspiration speed setting by pressing **ENTER**.
4. The down arrow appears on the display.
5. Use the ∇ keys to select dispensing speed.
6. Confirm settings by pressing **ENTER**. You are returned to the main menu.

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CHANGING MODE

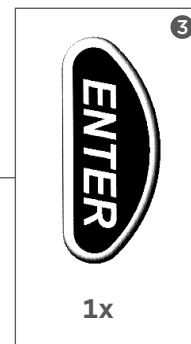
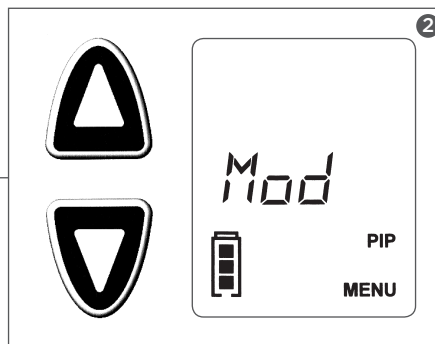
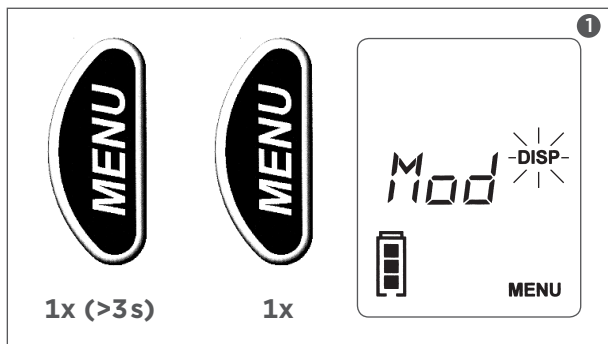
1. Hold down the **MENU** key (>3s) and then press it again briefly.
 - The mode can then be selected.
 - Make sure that plunger is at lowest position!
 - Modes available are **PIP** and **DISP**.
 - **DISP** is the default mode.
2. Use the $\nabla\Delta$ keys to select desired mode.
3. Confirm setting by pressing **ENTER**.

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USE OF COMPATIBLE TIPS

Compatible dispenser tips which are not manufactured by Gilson are not encoded for size. These tips can still be used with the REPETMAN (see **INSERTING TIPS**, on page 16). After insertion of one of these tips, the display will automatically go to tip selection mode.

By default, the 5 mL tip is displayed flashing regardless of the inserted tip. Ten possible tip sizes are available.

1. Use the $\nabla\Delta$ keys to select appropriate tip size.
2. Confirm setting by pressing **ENTER**.

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RESIDUAL STROKE

A lock prevents the residual stroke from being dispensed as a regular dispensed or pipetted volume. The symbol ▽ indicates this remaining volume. It can be discharged by holding the **STEP** key.

“FREE-JET DISPENSING”

When tips with volumes of 5 mL or larger are selected, the default dispensing speed adjusts to allow accurate dispensing without contacting the walls of vessel (free—jet dispensing). This can be overridden by selecting a different speed (page 28).

MAINTENANCE AND CLEANING

REPETMAN is calibrated at the factory and is maintenance-free. If required, clean the instrument with a cloth dampened with water or a mild soap solution. **REPETMAN must not be dismantled.**

DEFAULT SETTINGS

To revert REPETMAN to its factory default settings, simultaneously press the **MENU** and **ENTER** keys for >3 seconds, with the plunger in the lowest position.

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TROUBLESHOOTING

PROBLEM	ERROR MESSAGE DISPLAYED	POSSIBLE CAUSE	CORRECTIVE ACTION
Tip dripping	No message	Leaking tip	Replace tip
Faint display; reduced speed.	Battery symbol flashing	Battery low	Recharge battery pack or replace with charged battery pack.
Instrument not responding.	ERR 001	<ul style="list-style-type: none"> Locking lever not engaged during aspirating or dispensing. Locking lever open at lower third of the tip. 	<ol style="list-style-type: none"> Press ENTER to confirm error. <p>Follow MENU instructions to avoid instrument damage!</p> <p>Display: CLO (Close).</p> <ol style="list-style-type: none"> Engage the locking lever. <p>Display: rEF (Reference run).</p> <ol style="list-style-type: none"> Press the STEP key. <p>Attention: Liquid remaining in tip will be discharged during the reference run (rEF). Discharge sample into a suitable vessel (see pages 5, point 10).</p>
Instrument not responding.	ERR 002	<p>Motor blocked</p> <ul style="list-style-type: none"> Battery low. Viscosity is too high (see page 7). Tip orifice obstructed. 	<ol style="list-style-type: none"> Press ENTER to confirm error. <p>Follow MENU instructions to avoid instrument damage!</p> <ul style="list-style-type: none"> Recharge or replace battery pack. Reduce aspirating/dispensing speed. <p>Display: rEF (Reference run).</p> <ol style="list-style-type: none"> Press the STEP key <p>Attention: Liquid remaining in tip will be discharged during the reference run (rEF). Discharge sample into a suitable vessel (see pages 5, point 10).</p>
Instrument not responding.	ERR TIP	Locking lever opened too early within the lower third of the tip.	<ol style="list-style-type: none"> Press ENTER to confirm error. <p>Follow MENU instructions to avoid instrument damage!</p> <ol style="list-style-type: none"> Display: OPE (Open). Push the locking lever to its upper stop.

TROUBLESHOOTING - CONTINUED-

PROBLEM	ERROR MESSAGE DISPLAYED	POSSIBLE CAUSE	CORRECTIVE ACTION
Instrument not responding - continued -	-	-	<ol style="list-style-type: none"> 3. Display alternating between out TIP and CLO. Remove tip and close locking lever (without tip). 4. REPETMAN automatically carries out a reference run. 5. Display: no TIP Insert new tip (see page 16).
Instrument not responding.	ERR 003	Electronic fault or tip connector not at lowest position when inserting new tip.	<ol style="list-style-type: none"> 1. Press ENTER to confirm error. 2. Follow MENU instructions: <ul style="list-style-type: none"> • Display rEF (reference run). • Press STEP Key. <p>Attention: Liquid remaining in tip will be discharged during reference run (rEF). Discharge sample into suitable vessel (see page 5, point 10). If error message is still shown on the display remove battery pack and insert it again (reset). If this does not solve the problem send instrument back for repair.</p>
No tone, no light when placing the REPETMAN into the charging dock.	ACC	<ul style="list-style-type: none"> • Instrument contains no battery. • Battery pack inserted improperly. • Battery pack faulty. 	<ul style="list-style-type: none"> • Insert battery pack. • Make sure battery contacts are aligned with contacts. • Replace battery pack.
Charge—indicator is flashing	-	Removal or insertion of a new battery pack	<ul style="list-style-type: none"> • Charge battery pack for at least 3.5 hours inside the REPETMAN.
Dispensing process interrupted. Attention! Volume not completely discharged.	ERR VOL	Dispensing process interrupted by: <ul style="list-style-type: none"> • Pressing the Δ key during dispensing. • electronic faults. 	<ol style="list-style-type: none"> 1. Press ENTER to confirm error. 2. Restart dispensing by pressing the STEP key.

Note:

If other error messages are displayed, remove battery pack and insert it again. If this does not solve the problem, send instrument in for repair.

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TECHNICAL DATA

Battery pack:

Rechargeable Nickel Metal Hydride
battery pack:
NiMH, 4 x AAA cells,
4,8 V/700 mAh

Charging times for empty battery pack:

- in battery compartment of
charging dock: approx. 2.5h,
- inside the REPETMAN: approx.
3.5h,
- when battery capacity
starts to diminish, recharge
battery pack in the battery
compartment of the charging
dock.

Battery pack capacity:

A full charge gives approx. 2.5 hours
of continuous work with a 12.5 mL
REPET-TIP.

This sign certifies that the product
meets the requirements of the EC
directives (In Vitro Diagnostic
Medical Devices (98/79/EC), Safety
requirements for electrical
laboratory use, Low voltage
directive (2014/35/EU),
Electromagnetic Compatibility, EMC
(2014/30/EU)) and has been tested
according to the specified test
methods.

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RETURN FOR REPAIR

Caution!

For safety reasons, only clean and decontaminated instruments will be accepted for inspection or repair.

- Thoroughly clean and decontaminate the instrument.
- Send back the pipette in the safety bag delivered with it to your local distributor.

Return for repair:

Return shipment has to be effected at the risk and the cost of the sender.

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DISPOSAL

The adjoining symbol means that storage batteries and electronic devices must be disposed of separately from household trash (mixed municipal waste) at the end of their service life.

- According to the Directive 2002/96/EC of the European Parliament and of the Council on Waste Electrical and Electronic Equipment (WEEE) of 27 January 2003, electronic equipment requires disposal according to the relevant national disposal regulations.
- Batteries contain substances that can have harmful effects on the environment and human health. Therefore according to the Directive 2006/66/EC of the European Parliament and the Council on Waste Batteries of 6 September 2006 batteries require disposal according to the relevant national disposal regulations. Dispose of batteries only when completely discharged.

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Warning!

Do not short-circuit the battery to discharge it!

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WARRANTY

We shall not be liable for the consequences of improper handling, use, servicing, operation or unauthorized repairs of the instrument or the consequences of normal wear and tear especially of wearing parts such as pistons, seals, valves and the breakage of glass as well as the failure to follow the instructions of the operating manual. We are not liable for damage resulting from any actions not described in the operating manual or if non-original parts have been used.

Subject to technical modification without notice. Errors excepted.

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LT801491 | ©2017 Gilson, Inc. All rights reserved.